

# NEUATION

TECHNOLOGIES



## iSTIR MS10

### PRODUCT USER MANUAL

# TABLE OF CONTENT

1.	Introduction	1
2.	Intended Use	1
3.	Safety Instructions	1
4.	Technical Specification	2
5.	Features	3
6.	Accessories	3
7.	Standard Parts Listing	3
8.	Installation	4
9.	User Interface & Display	4
10.	Operation	5
11.	Troubleshooting	6
12.	Maintenance & Cleaning	7
13.	Warranty Statement	7
14.	Product Disposal	8

## 1. INTRODUCTION

This manual provides important safety information for this Magnetic Stirrer. It should be kept near the equipment for quick & easy reference. This stirrer is specially designed for precise control of stirring speed which allows gentle to vigorous mixing with speed ranging from 200 to 2200 RPM with a maximum capacity of 10 Ltr. The multipurpose digital display assists users by displaying various parameters like speed, time, pulse & mode of operation, etc.



## 2. INTENDED USE

The hotplate stirrer is suitable to use for mixing and/or heating the liquids with the maximum capacity of 10 Ltr. It is designed for use in general laboratories, pharmacies, schools and universities.

**NOTE:** Before using the instrument, please read this user manual carefully. This user manual is intended to assist with the operation and care of the unit only and not its repair. For repair please contact the supplier.

## 3. SAFETY INSTRUCTIONS

In order to use this equipment properly & safely, every user must read the user manual & observe safety instructions.

1. Place the device on a flat, stable, clean, non-slippery and fire-proof surface.
2. Ensure that only trained staff uses the device. Keep the instructions manual in a place where it can be accessed easily.
3. Beware of the possible effects of magnetic field on pacemakers, data media, etc.
4. Properly lift the device with both hands while moving or installing.
5. Do not place any steel or magnetic material on the top surface except  **WARNING** the recommended magnetic stirring bar with beaker or flask in between. Doing this can affect the magnetism of device.
6. Wear your personal protective equipment in accordance with the hazardous category of the media to be processed. Otherwise, there is a risk from:
  - ☞ Splashing of liquids.
  - ☞ Ejection of parts.
  - ☞ The release of toxic or combustible gases.
7. The top surface must be clean and kept intact. Wear protective gloves  **WARNING** while cleaning the device. Use a soft mild cloth to clean.
8. Do not use damaged beaker, flask, stirring bar or any other component for operation. It may affect the efficiency of the device.
9. Reduce the speed if:
  - ☞ If the medium splashes out of the vessel due to high speed.
  - ☞ If the appliance is not running smoothly.
  - ☞ If the container moves on the base plate.

### 3. SAFETY INSTRUCTIONS

10. Do not move the device when it's connected to power supply or during its operation.
11. Ensure that the product is used only for specified operation. It should not be used for shaking hazardous or reactive solutions.
12. Give specific attention to the risks associated with:
  - ☞ Flammable materials.
  - ☞ Flammable media with low steam pressure.
  - ☞ Overfilling of medium.
  - ☞ Incorrect vessel size.
  - ☞ Unstable vessel.
  - ☞ Glass breakage.
13. Note that, there is a possibility of contamination or unwanted chemical reaction.
14. The rating of power supply must match with the specified rating.
15. Process pathogenic materials only in a closed vessel under a suitable extractor hood. Do not operate the appliance in explosive atmospheres, with the hazardous substance or under water.
16. Only that liquid should be used for mixing which does not react dangerously to the extra energy produced through processing. This also applies to any extra energy produced in other ways. For example: Through light irradiation, through surrounding temperature etc.
17. Abrasion of the dispersion equipment or the rotating accessories can get into the medium you are working on.
18. The chemical reaction of PTFE can occur when in contact with molten or dissolved alkaline earth metals. Only elementary fluorine, chlorine trifluorine and alkaline metals do attack PTFE, halogen hydrocarbons have a reversible swelling effect.
19. To protect do not cover the device, even partially, with elements such as metallic plates or sheets otherwise it may overheat. Ensure that mounting plate remains clean.



### 4. TECHNICAL SPECIFICATIONS

Motor	DC
Max Stirring Quantity on Equip(H <sub>2</sub> O)	10 Liters
Speed range	200 to 2200 RPM
Speed control	10 RPM / Step
Accuracy	±20 RPM
Time range	1 to 999 minutes & Infinite
Bar Size	25-60 mm
Setup plate	SS



## 4. TECHNICAL SPECIFICATIONS

Dimensions (L x W x H)	156 x 248 x 77 mm
Protection class according to DIN EN	IP 43
Permissible ambient temp.	5 - 40°C
Permissible relative humidity	80%
Weight	1.4 Kg
Input Voltage	24V $\overline{\text{---}}$ 0.3A
Power Consumption	5.5 W

Maximum recommended bar size	Maximum Speed
25mm	2200 RPM
35mm	2200 RPM
50mm	1650 RPM
60mm	1500 RPM

## 5. FEATURES


- Powerful motor delivers constant speed under varying load conditions.
- Variable speed setting from 200 to 2200 RPM in steps of 10 RPM & long timer range from 1-999 mins & infinite mode.
- Microprocessor controlled with last run memory feature.
- Large & clear display shows critical parameters for easy reading.
- Small footprint saving valuable bench space.

## 6. ACCESSORIES


- Magnetic Stirring Bar.
- Power Adaptor: 100-240VAC 50/60Hz (Input),  
24V  $\overline{\text{---}}$  1.5A (Output).
- User Manual and Warranty Card.

## 7. STANDARD PARTS

Front Side

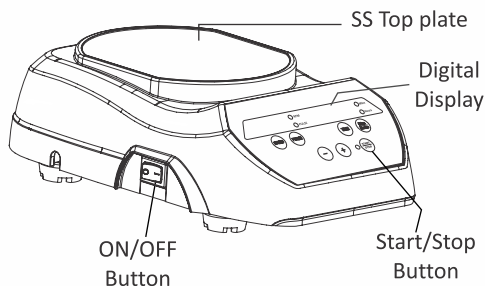

Rear Side

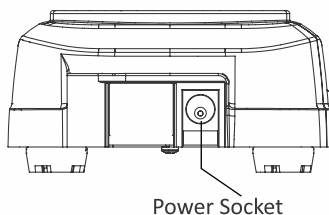

## 8. INSTALLATION

This stirrer is provided in a box. Open the box, then remove the packaging and gently place the device on the firm & leveled surface. Take care while unpacking & removing all accessories. The user manual should be kept with the device for easy access. Please keep all packaging in safe storage for at least two years for warranty purpose.


### 8.1 LOCATION & MOUNTING

Place the stirrer on a flat and leveled surface & ensure that all the four legs of this stirrer stands on the surface firmly. Avoid installing on a slippery surface or surface prone to vibration.


1. Ideal ambient temperature is  $25^{\circ}\text{C} \pm 1^{\circ}\text{C}$ ; avoid placing the unit in direct sunlight.
2. Keep clearance of at least 20 cm from all side to increase the cooling efficiency.
3. Keep the unit away from heated solution to avoid sample temperature issues.
4. Do not place the equipment at a place where it becomes difficult to operate it.

### 8.2 CONNECTING POWER ADAPTOR

1. Connect one end of the power adaptor to the rear side of the stirrer and another end to the power supply as shown in the figure below.
2. Push the power adaptor firmly for proper connection and turn ON the main switch.
3. Make sure that input power source is according to stirrer requirement.

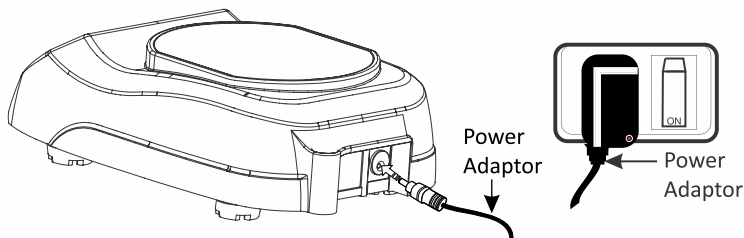

## 9. USER INTERFACE & DISPLAY

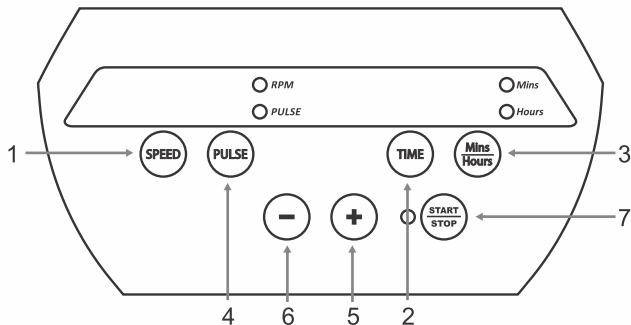

## 9. USER INTERFACE & DISPLAY


Display			
RPM	Mins	ACTUAL	Pulse
2200	999	200	30
Display shows current <b>speed</b> & <b>time</b> values		Display shows Speed, Time and Pulse Mode	

Item	Name	Function
1	Speed	Press "Speed" to select Speed value. Then press "+/-" to set Speed value in RPM
2	Time	Press "Time" to select time value. Then press "+/-" to set Time value in minutes
3	Mins/Hours	Press "Mins / Hours" to display time either in Minutes or Hour/s
4	Pulse	Press "Pulse" to activate the pulse mode in any operation. Long pressing "Pulse" lets user select timings for pulse mode from 30sec to 99sec.
5	Increment	Press "+" to increase values
6	Decrement	Press "-" to decrease values
7	Start/Stop	Press "Start/Stop" to Start or Stop any operation

## 10. OPERATION

### 10.1 SWITCHING ON THE DEVICE

Plug the power adaptor into its socket on the rear panel, now connect the power cord to AC power output. Now switch ON the mains and switch ON the stirrer from the front side, as shown in the figure. The power LED on its front panel will glow.

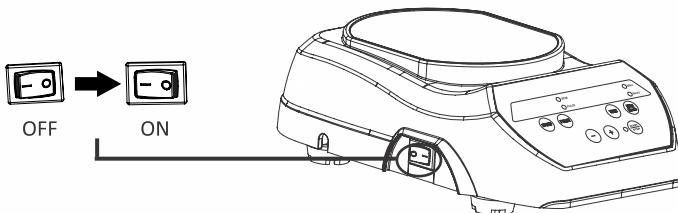

**Important Note:** Values of any parameter will get set saved only after the respective display blinks for 5 times.

## 10. OPERATION

### 10.2 SETTING TIME

The time in the display will show in minutes. The stirrer timer set for operation is in between 1min to 999 mins and infinite. Infinite time is seen as “∞”. To set time, press “TIME” button and then set the value by pressing “+/-” button. When time button is pressed, the “Mins” display will blink indicating the time value is selected. Once time value is selected, single press “+/-” will increase or decrease time value by 1min.

Long press increment or decrement button to quickly increase or decrease time value. Time value gets saved after multiple blinks. The timer is a countdown timer showing remaining time left. Once time is over, the stirrer stops.

### 10.3 SETTING SPEED

Speed is displayed as RPM. The minimum and maximum speed of stirrer is 200 RPM and 2200 RPM respectively. When the equipment is used for the 1<sup>st</sup> time, the speed will be set to zero and shown as “0”. Once speed value is selected, single press “+/-” will increase or decrease speed value by 10 RPM. Long press “+/-” to quickly increase or decrease speed value. Speed value gets saved after 5 blinks.

To operate the device as a heater set the time and temperature as informed in the previous section and set the speed of the stirrer to zero by pressing “+/-”.

### 10.4 PULSE MODE OPERATION

User can also program the pulse mode and set it for alternating times from 30 secs to 99 secs.

First set the required speed and total run time for operation. Then hold down the “Pulse” button to select the cycle time of stirring. The cycle time value will appear on the right side display, “SEC” will appear in left side display and pulse LED will start blinking indicating that the user can now modify the cycle time. Use increment or decrement button to modify the cycle time between 30 seconds to 99 seconds. Once the cycle time is modified, it will automatically save after 5 blinks. Press the “Pulse Button” once to START pulse mode operation with the modified cycle time.

Press “Pulse” again to stop pulse mode operation. User can also Press “Start / Stop” to STOP ongoing pulse mode operation.

## 11. TROUBLESHOOTING

**Any type of malfunction during operation can be identified by error an message on display. Proceed as follows in such cases.**

### 1. Problem : No Display.

**Root Cause : No main power connection.**

**Solution :** 1) Check main power is available.

## 11. TROUBLESHOOTING

- 2) Check that power adaptor is working properly and it is properly connected at both ends.
- 3) If the problem still persists then call authorized service person.

### 2. Problem : Err 2.

**Root Cause : Motor blockage/over load/fail, Encoder sensor fail.**

Solution : 1) First switch off the equipment, plug it out and Plug it in again.

- 2) If the problem still persists then call authorized service person.

## 12. MAINTENANCE & CLEANING

1. Before cleaning the device, allow the instrument to attain room temperature remove the power cord from the mains.
2. To clean the housing of the device from dyes, building materials or cosmetic materials use isopropyl alcohol as a cleaning agent.
3. To clean the housing of the device from food materials or fuels, use water containing detergent as cleaning agent.
4. Wear proper protective gloves while cleaning the instrument.
5. Be careful that no liquid enters the device during cleaning.
6. Kindly contact suppliers before using any other methods for cleaning.

## 13. WARRANTY STATEMENT

This product is warranted to be free from defects in material and workmanship for a period of two (2) years from the date of purchase. Your product will be duly repaired upon prompt notification in compliance with the following conditions :

This warranty is valid only if the product is used for its intended purpose and within the guidelines specified in this instruction manual. This warranty does not cover damage caused by accident, neglect, misuse, improper service, natural forces or other causes not arising from defects in original material or workmanship. This warranty does not cover any incidental or consequential damages, commercial loss or any other damages from the use of this product.

The warranty is invalidated by any non-factory modification, which will immediately terminate all liabilities on us for the products or damages caused by its use. The buyer and its customer shall be responsible for the product or use of products as well as any supervision required for safety. If requested the products must be returned to the distributor in well packed and insured manner and all shipping charges must be paid.

Some states do not allow limitation on the length of implied warranties or the exclusion or limitation of incidental or consequential damages. This warranty gives you specific legal rights. This warranty is given expressly in lieu of all other warranties, expressed or implied.

### 13. WARRANTY STATEMENT

The purchaser agrees that there is no warranty of merchantability or of fitness for any intended purpose and that there are no other remedies or warranties, expressed or implied, which extend beyond the description on the face of the agreement. This warranty is only applicable to the original purchaser.

Products received without proper authorization will not be entertained. All items returned for service should be sent postage prepaid in the original packaging or other suitable carton, padded to avoid damage. We will not be responsible for damage incurred by improper packaging.

All items returned for service should be set postage prepaid in the original packaging or other suitable carton, added to avoid damage.

This warranty is valid only if the warranty is registered with the supplier within 30 days from the date of purchase.

For you reference, make a note of serial number, date of purchase and supplier here.	
Serial No.:	Purchase Date:
Supplier:	

### 14. PRODUCT DISPOSAL



In case the product is to be disposed of, the relevant legal regulations are to be observed.

Information on the disposal of electrical and electronic devices in the European Community

The disposal of electrical devices is regulated within the European Community by national regulations based on EU Directive 2012/19/EU on waste electrical and electronic equipment (WEEE). According to these regulations, any devices supplied after 13.06.05 in the business to the business sphere, to which this product is assigned, may no longer be disposed of in municipal or domestic waste. They are marked with the following symbol to indicate this.

As disposal regulations within the EU may vary from country to country, please contact your supplier if necessary.

## Neuation Technologies Pvt. Ltd.

Plot No. 15, GIDC Electronic Park SEZ Kolavada Road, Gandhinagar – 382026, Gujarat, India

---

**Website:** [www.neuation.com](http://www.neuation.com)